

Manual

EpiScript RNase H– Reverse Transcriptase

For Laboratory Use.

Manual

EpiScript RNase H- Reverse Transcriptase

Contents

1. Intended use	3
2. Warning and limitations	3
3. Product designations	4
4. Product specifications	4
5. General considerations for cDNA synthesis	4
6. EpiScript RNase H- Reverse Transcriptase cDNA synthesis procedure	6
7. Further support	7

Manual

EpiScript RNase H– Reverse Transcriptase

1. Intended use

EpiScript™ RNase H– Reverse Transcriptase (EpiScript RT) is a recombinant MMLV reverse transcriptase with reduced RNase H activity. It is active up to 55 °C to achieve higher specificity and is efficient at producing full-length cDNA from long RNA templates.

EpiScript RT demonstrates significantly more activity than other MMLV Reverse transcriptase enzymes and is capable of producing cDNA from as little as 50 pg of total RNA for real-time RT-PCR (qRT-PCR) analysis and other applications.

EpiScript RT is intended for use as a general purpose reagent in molecular diagnostic assays that are based on RT-PCR target detection technologies by clinical laboratory professionals located in the United States. This product must be qualified and validated by clinical laboratory end users for suitability in the detection of any specific target using the procedures in this Manual as guidance.

EpiScript RT is supplied at a concentration of 200 U/μL.

2. Warning and limitations

This product, when used in combination with nucleic acid amplification master mixes and oligonucleotide primers and probes for specific targets, may be used as a component in molecular diagnostic assays upon qualification and validation by the clinical laboratory end-user.

This product is classified as a general purpose *in vitro* diagnostic device reagent in the United States as defined in: <https://www.accessdata.fda.gov/scripts/cdrh/cfdocs/cfcfr/CFRSearch.cfm?FR=864.4010>



Medical Device Manufacturer
LGC Biosearch Technologies
2905 Parmenter St.
Middleton, WI USA 53562

Manual

EpiScript RNase H- Reverse Transcriptase

3. Product designations

Product	Size	Catalogue number	Reagent description	Part number	Volume
EpiScript RNase H- Reverse Transcriptase Kit	10,000 units	ERT12910K	EpiScript RNase H- Reverse Transcriptase, 200 U/ μ L	E0144-200D1	50 μ L
			10X RT Reaction Buffer	SS000737-D2	250 μ L
			DTT (100 mM)	SS000065-D6	250 μ L
EpiScript RNase H- Reverse Transcriptase Kit	25,000 units	ERT12925K	EpiScript RNase H- Reverse Transcriptase, 200 U/ μ L	E0144-200D2	125 μ L
			10X RT Reaction Buffer	SS000737-D3	600 μ L
			DTT (100 mM)	SS000065-D8	600 μ L

Table 1. EpiScript RNase H- Reverse Transcriptase Kit products and codes.

Product	Size	Catalogue number	Reagent description	Part number	Volume
EpiScript RNase H- Reverse Transcriptase	25,000 units	ERT12925K-ENZ	EpiScript RNase H- Reverse Transcriptase, 200 U/ μ L	E0144-200D2	0.125 mL
	250,000 units	ERT12925K-1.25mL	EpiScript RNase H- Reverse Transcriptase, 200 U/ μ L	E0144-200D3	1.25 mL
	2,500,000 units	ERT12925K-12.5mL	EpiScript RNase H- Reverse Transcriptase, 200 U/ μ L	E0144-200D4	12.5 mL

Table 2. EpiScript RNase H- Reverse Transcriptase standalone enzyme products and codes.

4. Product specifications

Storage: Store only at -20 °C in a freezer without a defrost cycle.

Storage buffer: EpiScript RT is supplied in a 50% glycerol solution containing 50 mM Tris- HCl (pH 7.5), 100 mM sodium chloride, 1 mM DTT, 0.1 mM EDTA and 0.1% Triton® X-100.

Unit definition: One unit of EpiScript RT catalyses the incorporation of 1 nmol of dTTP into acid-insoluble material in 10 minutes at 37 °C using saturating amounts of oligo(dT)-primed poly(rA) as template.

Contaminating activity assays: All EpiScript RT components are free of detectable contaminating DNA exonuclease and endonuclease, and RNase activities.

5. General considerations for cDNA synthesis

Maintaining an RNase-free environment:

Ribonuclease contamination is a significant concern for those working with RNA. The ubiquitous RNase A is a highly stable and active ribonuclease that can contaminate any lab environment and is present on human skin. However, creating an RNase-free work environment and maintaining RNase-free solutions is critical for performing successful cDNA synthesis reactions. Therefore, we strongly recommend that the user:

Manual

EpiScript RNase H- Reverse Transcriptase

- 1) Autoclave all tubes and pipette tips that will be used in the cDNA synthesis reactions.
- 2) Always wear gloves when handling samples containing RNA. Change gloves frequently especially after touching potential sources of RNase contamination such as door knobs, pens, pencils and human skin.
- 3) Always wear gloves when handling kit components. Do not pick up any kit component with an ungloved hand.
- 4) Keep all kit components tightly sealed when not in use. Keep all tubes containing RNA tightly sealed during the incubation steps.

We *strongly* recommend the addition of LGC Biosearch Technologies' RiboGuard™ RNase Inhibitor to each reaction.

Choice of primer for first-strand cDNA synthesis:

First-strand cDNA synthesis can be primed using either an **oligo(dT) primer**, **random primers**, or **gene-specific primers** (all primers to be provided by the user).

An **oligo(dT) primer** is the most commonly used method for priming first-strand cDNA synthesis when using an eukaryotic RNA sample. Oligo(dT) primes cDNA synthesis only from the poly(A) tail present at the 3' end of almost all eukaryotic mRNAs. Since poly(A) RNA constitutes just 1-5% of the RNA in a eukaryotic total cellular RNA preparation, the complexity of the cDNA produced is significantly less than when the cDNA is synthesised using random primers. Lower-complexity cDNA can result in a more sensitive and specific PCR amplification. Additionally, priming cDNA synthesis with an oligo(dT) primer precludes the need to enrich the RNA sample for poly(A) RNA. Typically, the oligo(dT) primer is 18-21 nucleotides in length.

Random primers initiate ccDNA synthesis from all RNA species (rRNA and mRNA) in a total cellular RNA sample. Since rRNA, which constitutes >95% of the RNA in a total RNA sample, is converted to cDNA using random primers, the complexity of the cDNA will be much greater than when priming the reaction with oligo(dT). As a result, random primers are much less frequently used than oligo(dT) primers. Random primers are typically 6-9 nucleotides in length and can be helpful when:

- 1) Synthesising cDNA from mRNAs that lack a poly(A) tail (such as bacterial mRNA) or have a very short poly(A) tail;
- 2) Priming cDNA synthesis from partially degraded RNA samples such as those obtained from formalin-fixed paraffin-embedded (FFPE) tissue sample;
- 3) Priming cDNA synthesis of a poly(A)-enriched RNA sample;
- 4) It is necessary to eliminate or reduce 3' sequence bias that can result when using an oligo(dT) primer.

Gene-specific primers, designed and synthesised by the user, provide the greatest specificity when priming cDNA synthesis of an mRNA. However, the user frequently must determine the optimal primer annealing and extension (reverse transcription) conditions empirically for each primer used.

Manual

EpiScript RNase H- Reverse Transcriptase

6. EpiScript RNase H- Reverse Transcriptase cDNA synthesis procedure

The following protocol has been optimised to convert 100 pg of cellular RNA to 1 µg of first-strand cDNA using an oligo(dT) primer or random primers. The use of gene-specific primers may require additional optimisation of the reaction.

Gently mix and briefly centrifuge all kit components prior to dispensing.

1. Denature the RNA sample and anneal the primer(s). For each reaction, combine the following components on ice, in a sterile (RNase-free) 0.2 mL or 0.5 mL tube:

x	µL RNase-free water
x	µL total RNA sample (up to 1 µg)
10	pmol oligo(dT) ₁₈₋₂₁ primer
	- or -
100	ng random primers
	- or -
x	µL gene-specific primers
<hr/>	
10	µL total reaction volume

2. Incubate at 65 °C for 2 minutes in a water bath or thermal cycler with heated lid.
3. Chill on ice for 1 minute. Centrifuge briefly in a microcentrifuge.
4. For each reaction, combine the following components on ice:

x	µL RNase-free water	
10	µL annealed RNA:primer(s) (from step 3 above)	
2	µL 10X RT Reaction Buffer	
2	µL 100 mM DTT	
0.5	µL RiboGuard RNase Inhibitor (optional, sold separately)	
1	µL 10 mM dATP	Solution (provided by the user, sold separately)
1	µL 10 mM dCTP	Solution (provided by the user, sold separately)
1	µL 10 mM dGTP	Solution (provided by the user, sold separately)
1	µL 10 mM dTTP	Solution (provided by the user, sold separately)
0.5	µL EpiScript Reverse Transcriptase (200 U/µL)	
<hr/>		
20	µL total reaction volume	

Manual

EpiScript RNase H- Reverse Transcriptase

5. Mix the reaction gently.
If using oligo(dT) primers, incubate the reaction at 37 °C for 60 minutes.
If using random primers, incubate the reaction at room temperature for 10 minutes and then at 37 °C for 60 minutes.
6. Terminate the reaction by heating at 85 °C for 5 minutes.
7. Chill on ice for at least 1 minute. Centrifuge briefly in a microcentrifuge.
8. The cDNA can be used immediately, without purification, for end-point or real-time PCR (qPCR), converted to double-stranded cDNA, or stored at -20 °C for future use.

7. Further support

If you require any further support, please do not hesitate to contact our Technical Support Team: techsupport@lgcgroup.com.



For Laboratory Use.

Integrated tools. Accelerated science.

   @LGCBiosearch | biosearchtech.com

All trademarks and registered trademarks mentioned herein are the property of their respective owners. All other trademarks and registered trademarks are the property of LGC and its subsidiaries. Specifications, terms and pricing are subject to change. Not all products are available in all countries. Please consult your local sales representative for details. No part of this publication may be reproduced or transmitted in any form or by any means, electronic or mechanical, including photocopying, recording or any retrieval system, without the written permission of the copyright holder. © LGC Limited, 2022. All rights reserved. GEN/784/EK/0722



BIOSEARCH™
TECHNOLOGIES