

# Manual

## T4 Polynucleotide Kinase, Cloned

For Research Use Only. Not for use in diagnostic procedures.

T4 Polynucleotide Kinase, Cloned is part of the Epicentre™ product line, known for its unique genomics kits, enzymes, and reagents which offer high quality and reliable performance.

**BIOSEARCH**  
TECHNOLOGIES™  
GENOMIC ANALYSIS BY LGC

# Manual

T4 Polynucleotide Kinase, Cloned

---

## Contents

<a href="#"><u>1. Introduction</u></a> .....	<a href="#"><u>3</u></a>
<a href="#"><u>2. Product designations and kit components</u></a> .....	<a href="#"><u>3</u></a>
<a href="#"><u>3. Product specifications</u></a> .....	<a href="#"><u>3</u></a>
<a href="#"><u>4. Protocol I: end-labeling with T4 PNK</u></a> .....	<a href="#"><u>4</u></a>
<a href="#"><u>5. Protocol II: phosphorylation of oligodeoxynucleotides with T4 PNK</u></a> .....	<a href="#"><u>4</u></a>
<a href="#"><u>6. Further support</u></a> .....	<a href="#"><u>5</u></a>

# Manual

## T4 Polynucleotide Kinase, Cloned

### 1. Introduction

T4 Polynucleotide Kinase, Cloned (T4 PNK) catalyses the transfer of the  $\gamma$ -phosphate of ATP to the 5' terminus of single- and double-stranded DNA or RNA molecules that have a 5' hydroxyl. The enzyme also removes the 3' phosphate from 3'-phosphoryl polynucleotides, deoxyribonucleoside 3'-monophosphates and deoxyribonucleoside 3',5'-diphosphates to form a 3' hydroxyl. T4 PNK has many uses, including labeling the 5' ends of DNA or RNA with  $^{32}\text{P}$  for use in sequencing, as hybridisation probes or in transcript mapping with S1 nuclease or Mung Bean Nuclease. The enzyme can also be used to phosphorylate oligodeoxynucleotide linkers or other DNA molecules prior to ligation and to end-label oligodeoxynucleotide primers for use in sequencing reactions. T4 PNK may also be used in an exchange reaction to radiolabel the 5' ends of DNA or RNA molecules that have unlabeled 5' phosphates. It also may be used to dephosphorylate the 3'-ends of RNA in the absence of ATP.

### 2. Product designations and kit components

Product	Kit size	Catalogue number	Reagent description	Part number	Volume
T4 Polynucleotide Kinase, Cloned	3,000 units	P0503K	T4 PNK (10 U/ $\mu\text{L}$ )	E0055-10D	300 $\mu\text{L}$
			10X T4 PNK Buffer	SS000272-D7	1.5 mL

### 3. Product specifications

**Storage:** Store only at -20 °C in a freezer without a defrost cycle.

**Storage buffer:** T4 PNK is supplied in a 50% glycerol solution containing 50 mM Tris-HCl (pH 7.5), 0.1 M NaCl, 0.1 mM EDTA, 1 mM dithiothreitol (DTT) and 0.1% Triton® X-100 (Rohm & Haas).

**Unit definition:** One unit catalyses the transfer of 1 nmol of phosphate from  $\gamma$ -[ $^{32}\text{P}$ ]-labeled ATP to the 5'-hydroxyl ends of micrococcal nuclease-treated calf thymus DNA in 30 minutes at 37 °C.

**10X T4 PNK Buffer:** 330 mM Tris-acetate (pH 7.5), 660 mM potassium acetate, 100 mM magnesium acetate and 5 mM DTT.

**10X T4 PNK Buffer does not contain ATP. A 100 mM ATP Solution may be purchased separately (LGC, Biosearch Technologies™ RA02825).**

**Quality control:** T4 PNK is function-tested in a reaction containing 50 mM Tris-HCl (pH 7.5), 10 mM MgCl<sub>2</sub>, 10 mM  $\beta$ -mercaptoethanol, 0.1 mM  $\gamma$ -[ $^{32}\text{P}$ ]-ATP, 0.2 mg/mL of micrococcal nuclease-treated calf thymus DNA and varying amounts of enzyme.

**Contaminating activity assays:** T4 PNK is free of detectable DNA exo- and endonuclease, and RNase activities.

# Manual

## T4 Polynucleotide Kinase, Cloned

### 4. Protocol I: end-labeling with T4 PNK

The following protocol was developed for end-labeling primers for use in many applications requiring end-labeled DNA.

1. Combine the following reaction components on ice in the order given:

	<u>Final concentration</u>
x µL deionised water	---
12.5 µL pmol primer	0.5 pmol/µL
2.5 µL 10X T4 PNK Reaction Buffer	1X
0.5 µL γ-[ <sup>32</sup> P]-ATP (80 µCi; 12 pmol; > 7,000 Ci/mmol)	~0.5 pmol/µL
3 U T4 PNK	0.12 U/µL
25 µL Total reaction volume	

2. Incubate at 37 °C for 30 minutes.
3. Inactivate the T4 PNK by incubating at 70 °C for 5 minutes.

End-labelled primers can be used in cycle sequencing reactions without separating unincorporated nucleotides from the labeled primer. Primers should be stored at -20 °C and may be used for up to 2 weeks after labelling.

### 5. Protocol II: phosphorylation of oligodeoxynucleotides with T4 PNK

The following protocol describes the nonradioactive phosphorylation of oligodeoxy- nucleotides used in ligation amplification assays with Ampligase™ Thermostable DNA Ligase. The protocol can also be used for other applications requiring the addition of a phosphate to the 5' end of a DNA strand.

1. Combine the following reaction components on ice in the order given:

	<u>Final concentration</u>
x µL deionised water	---
10 µL oligonucleotide at 1.0 µg/µL*	0.2 µg/µL
5 µL 10X T4 PNK Reaction Buffer	1X
5 µL 10 mM ATP Solution	1 mM
3 U T4 PNK	0.06 U/µL
50 µL Total reaction volume	

# Manual

## T4 Polynucleotide Kinase, Cloned

---

2. Incubate at 37 °C for 30 minutes.
3. Inactivate the T4 PNK by incubating at 70 °C for 5 minutes.

\*To determine the micrograms of double-stranded DNA needed in other applications, multiply the length of the DNA fragment by 0.00034. For example, add 1.7 µg of a 5,000 bp restriction fragment ( $5,000 \times 0.00034 = 1.7$ ).

### **6. Further support**

If you require any further support, please do not hesitate to contact our Technical Support Team:  
[techsupport@lgcgroup.com](mailto:techsupport@lgcgroup.com).

# Integrated tools. Accelerated science.

 @LGCBiosearch | [biosearchtech.com](http://biosearchtech.com)

All trademarks and registered trademarks mentioned herein are the property of their respective owners. All other trademarks and registered trademarks are the property of LGC and its subsidiaries. Specifications, terms and pricing are subject to change. Not all products are available in all countries. Please consult your local sales representative for details. No part of this publication may be reproduced or transmitted in any form or by any means, electronic or mechanical, including photocopying, recording or any retrieval system, without the written permission of the copyright holder. © LGC Limited, 2021. All rights reserved. GEN/884/EK/0821

For Research Use Only. Not for use in diagnostic procedures.



**BIOSEARCH**  
**TECHNOLOGIES**  
GENOMIC ANALYSIS BY LGC