

# Manual

## TransforMax EC100D *pir+* Electrocompetent *E. coli* and TransforMax EC100D *pir-116* Electrocompetent *E. coli*

---

For Research Use Only. Not for use in diagnostic procedures.

TransforMax™ EC100D™ *pir+* Electrocompetent *E. coli* TransforMax™ EC100D™ *pir-116* and Electrocompetent *E. coli* are part of the Epicentre™ product line, known for its unique genomics kits, enzymes, and reagents which offer high quality and reliable performance.

# Manual

TransforMax EC100D *pir+* and *pir-*116 Electrocompetent *E. coli*

---

## Contents

<a href="#">1. Product description</a> .....	3
<a href="#">2. Product specifications</a> .....	3
<a href="#">3. Product designations and kit components</a> .....	4
<a href="#">4. Storage conditions</a> .....	4
<a href="#">5. Electroporation procedure</a> .....	4
<a href="#">6. References</a> .....	5
<a href="#">7. Technical support and product guarantee</a> .....	5

# Manual

## TransforMax EC100D *pir+* and *pir*-116 Electrocompetent *E. coli*

### 1. Product description

The TransforMax EC100D *pir+* and the TransforMax EC100D *pir*-116 Electrocompetent *E. coli* each express the  $\Pi$  protein (*pir* gene product) for replication of vectors containing the R6K $\gamma$  conditional origin of replication (R6K $\gamma$ ori). The cells are derived from LGC, Biosearch Technologies' TransforMax EC100™ *E. coli* by P1 phage transduction with a strain containing either the *pir* or *pir*-116 gene linked to a dihydrofolate reductase (DHFR) marker. Both types of cells can be used for 'rescue cloning' of genomic DNA fragments transposed with an EZ-Tn5™ <R6K $\gamma$ ori/KAN-2> Transposon using either the EZ-Tn5 <R6K $\gamma$ ori/KAN-2> Insertion Kit or the EZ-Tn5 <R6K $\gamma$ ori/KAN-2> Tnp Transposome Kit. The TransforMax EC100D *pir+* Electrocompetent *E. coli* will maintain R6K $\gamma$ ori-containing vectors at approximately 15 copies per cell<sup>1</sup> for cloning of potentially toxic or unstable DNA sequences. The TransforMax EC100D *pir*-116 Electrocompetent *E. coli* are for high copy number propagation of up to 250 rescue plasmid copies per cell.<sup>1</sup>

### Important phenotypes and applications

- Expresses the  $\Pi$  protein (*pir* gene product) for rescue cloning and propagation of vectors containing an R6K $\gamma$  origin of replication.
- Compatible with vectors expressing the *LacZ'*  $\alpha$ -complementing peptide for 'blue/white' screening of recombinants.
- Restriction minus for efficient cloning of methylated genomic DNA.
- Accepts large clones for unbiased propagation and stability of large rescue clones.
- Endonuclease minus (*endA1*) to ensure high yields of plasmid clones.

### 2. Product specifications

	TransforMax EC100D <i>pir+</i> Electrocompetent <i>E. coli</i>	TransforMax EC100D <i>pir</i> -116 Electrocompetent <i>E. coli</i>
<b>Transformation efficiency:</b>	Greater than $5 \times 10^9$ cfu/ $\mu$ g of supercoiled DNA.	Greater than $5 \times 10^9$ cfu/ $\mu$ g of supercoiled DNA.
<b>Genotype</b>	F- <i>mcrA</i> $\Delta$ ( <i>mrr</i> - <i>hsdRMS</i> - <i>mcrBC</i> ) $\phi$ 80d <i>lacZ</i> $\Delta$ M15 $\Delta$ <i>lacX74</i> <i>recA1</i> <i>endA1</i> <i>araD139</i> $\Delta$ ( <i>ara</i> , <i>leu</i> )7697 <i>galU</i> <i>galK</i> $\lambda$ - <i>rpsL</i> <i>nupG</i> <i>pir</i> *(DHFR) Maintains plasmids at ~15 copies per cell.	F- <i>mcrA</i> $\Delta$ ( <i>mrr</i> - <i>hsdRMS</i> - <i>mcrBC</i> ) $\phi$ 80d <i>lacZ</i> $\Delta$ M15 $\Delta$ <i>lacX74</i> <i>recA1</i> <i>endA1</i> <i>araD139</i> $\Delta$ ( <i>ara</i> , <i>leu</i> )7697 <i>galU</i> <i>galK</i> $\lambda$ - <i>rpsL</i> <i>nupG</i> <i>pir</i> -116(DHFR) Maintains plasmids at ~250 copies per cell.
<b>Quality control:</b>	TransforMax EC100D <i>pir+</i> and <i>pir</i> -116 Electrocompetent <i>E. coli</i> yield $>5 \times 10^9$ transformants/ $\mu$ g of supercoiled DNA. Transformation efficiency is determined using 10 pg of pR6Kan control DNA, an Eppendorf Multipipettor (2.5 kV, fast charge rate), and a 2-mm cuvette. TransforMax EC100D <i>pir+</i> and <i>pir</i> -116 Electrocompetent <i>E. coli</i> are tested to free of contaminating DNA rendering resistance to ampicillin, tetracycline, kanamycin and chloramphenicol.	

# Manual

TransforMax EC100D *pir+* and *pir-116* Electrocompetent *E. coli*

## 3. Product designations and kit components

Product	Kit size	Catalogue number	Reagent description	Part number	Volume
TransforMax EC100D <i>pir+</i> Electrocompetent <i>E. coli</i>	5 x 100 µL*	ECP09500	TransforMax EC100D <i>pir+</i> Electrocompetent <i>E. coli</i>	SS001014-D	5 x 100 µL
			pR6Kan Control Plasmid**	SS000497-D	10 µL
TransforMax EC100D <i>pir-116+</i> Electrocompetent <i>E. coli</i>	5 x 100 µL*	EC6P095H	TransforMax EC100D <i>pir-116</i> Electrocompetent <i>E. coli</i>	SS001013-D	5 x 100 µL
			pR6Kan Control Plasmid**	SS000497-D	10 µL

\* 10 Electroporations.

\*\* Each is supplied with 10 µL (100 pg/µL) of pR6Kan Control DNA in TE Buffer.

## 4. Storage conditions

Store TransforMax EC100D *pir+* and *pir-116* *E. coli* cells at -70 °C and the pR6Kan Control DNA at either -20 °C or -70 °C.

## 5. Electroporation procedure

**NOTE:** The electroporation procedure described here uses 50 µL of electrocompetent cells. A different volume of cells can also be used based on the experiences and needs of the user. It is critical that the DNA, cuvettes, electroporation chamber (if applicable) and microcentrifuge tubes are thoroughly chilled on ice.

**CAUTION:** DNA should be prepared in water or very low ionic strength buffer (e.g. TE Buffer: 10 mM Tris-HCl [pH 7.5], 1 mM EDTA) to prevent arcing during electroporation. DNA from a Biosearch Technologies Fast-Link DNA ligation reaction can be diluted and used directly, without desalting or ethanol precipitation.

1. Thaw TransforMax EC100D *pir+* or *pir-116* cells on ice. Mix gently. Use the cells immediately. Unused cells can be refrozen at -70 °C.  
**NOTE:** Refrozen cells may have reduced transformation efficiency, but for most cases, the reduction is not significant enough to interfere with the desired results.
2. Transfer the desired amount of DNA and 50 µL of cells to a pre-chilled microcentrifuge tube.  
**NOTE:** A smaller volume of cells can be used based on the needs and experiences of the user. Mix the cells and DNA by pipetting up and down 2-3 times.  
**Control (optional):** Dilute the pR6Kan control DNA 1:10 with sterile, distilled water. Add 1 µL (10 pg) to 50 µL of cells as described.
3. Transfer the cell/DNA mix to the electroporation cuvette. Be sure that there are no air bubbles in the cuvette. Wipe the cuvette of any condensation. Place into the electroporator and apply the electric pulse at the manufacturer's recommendations for bacterial (*E. coli*) electroporation.
4. Immediately after electroporation, add 950 µL of room temperature SOC medium<sup>2</sup> to the cuvette. Mix gently by pipetting up and down 2-3 times.

# Manual

TransforMax EC100D *pir+* and *pir-*116 Electrocompetent *E. coli*

---

5. Transfer the cells to a sterile culture tube and incubate at 37 °C with shaking at 220-230 rpm for one hour to recover the cells and allow expression of the antibiotic resistance marker.
6. We recommend diluting the cells 1:10 and 1:100 and plating 100 µL of each dilution to the appropriate medium and antibiotic plate. Incubate the plates at 37 °C overnight.

**Control (optional):** Dilute the control reaction 1:20 and plate 100 µL (equivalent to 0.05 pg DNA) to LB-kanamycin (50 µg/mL) plates. If 250 colonies are observed on the plate, the transformation efficiency is  $5 \times 10^9$  cfu/µg or  $[(250 \text{ cfu}/0.05 \text{ pg DNA}) \times (10^6 \text{ pg}/\mu\text{g})]$ .

## 6. References

1. Metcalf WW *et al.*, (1994) *Gene* 138, 1.
2. Hanahan D, 1983, *J. Mol. Biol.*, 166, 557.

## 7. Further support

If you require any further support, please do not hesitate to contact our Technical Support Team: [techsupport@lgcgroup.com](mailto:techsupport@lgcgroup.com).



# Integrated tools. Accelerated science.

   @LGCBiosearch | [biosearchtech.com](https://biosearchtech.com)

*All trademarks and registered trademarks mentioned herein are the property of their respective owners. All other trademarks and registered trademarks are the property of LGC and its subsidiaries. Specifications, terms and pricing are subject to change. Not all products are available in all countries. Please consult your local sales representative for details. No part of this publication may be reproduced or transmitted in any form or by any means, electronic or mechanical, including photocopying, recording or any retrieval system, without the written permission of the copyright holder. © LGC Limited, 2021. All rights reserved. GEN/1016/EK/1121*

For Research Use Only. Not for use in diagnostic procedures.



**BIOSEARCH™**  
**TECHNOLOGIES**

GENOMIC ANALYSIS BY LGC