### **Product insert**



## **NxSeq Single-cell RNA-seq Beads**

For Research Use Only. Not for use in diagnostic procedures.

#### 1. Introduction

The NxSeq™ Single-cell RNA-seq Beads contain a covalently attached library of barcoded oligo-dT capture oligonucleotides to enable single-cell mRNA sequencing experiments. Every bead-attached oligo bead contains a 12 base, bead-specific barcode, a 14 base unique molecular identifier (UMI) that is distinct between oligos, and a 3' 30 base oligo-dT mRNA capture sequence. The structure of each bead and oligo sequence is as follows:

Bead - Linker -

J represents the bead barcode

N represents the 13 base UMI

V represents a non-dT base at position 14 of the UMI

T<sub>30</sub> represents the 30 base oligo-dT capture sequence

Please note that the NxSeq Beads contain an extended UMI of N13V (14 bases) compared to other commercially available beads. As such, adjust Read 1 length of your RNA sequencing accordingly.

#### 2. Product designations and kit components

Cat no.	Description	Size	Unit
NX-SCB-1	NxSeq Single-cell RNA-seq Beads	33 mg, 3 million beads	Each

#### 3. Bead re-suspension

When re-suspending the beads, use gentle flicking or pipetting of the solution. DO NOT VORTEX. Prior to use, wash the NxSeq Beads twice with 1 mL 100% ethanol, and then twice with 1 mL TE/TW (10 mM Tris pH 8.0, 1 mM EDTA, 0.01% Tween). Re-suspend the washed beads to the desired concentration in TE/TW.

#### 4. Recommended storage conditions

Store dry beads at -15 to -25 °C. Store beads re-suspended in TE/TW at +2 to +8 °C. Re-suspended beads are stable for one month.

# Integrated tools. Accelerated science.



lucigen.com biosearchtech.com

All trademarks and registered trademarks mentioned herein are the property of their respective owners. All other trademarks and registered trademarks are the property of LGC and its subsidiaries. Specifications, terms and pricing are subject to change. Not all products are available in all countries. Please consult your local sales representative for details. No part of this publication may be reproduced or transmitted in any form or by any means, electronic or mechanical, including photocopying, recording or any retrieval system, without the written permission of the copyright holder. © LGC Limited, 2020. All rights reserved. GEN/775/SW/1120

